


Serum and Dried Blood Spot (DBS) samples were compared using 33 matched samples in the range of 0 – 25 ng/mL. Passing Bablok analysis of the results yielded the following regression
 DBS AMH = 0 + 1.072 Serum AMH
 (r=0.975)

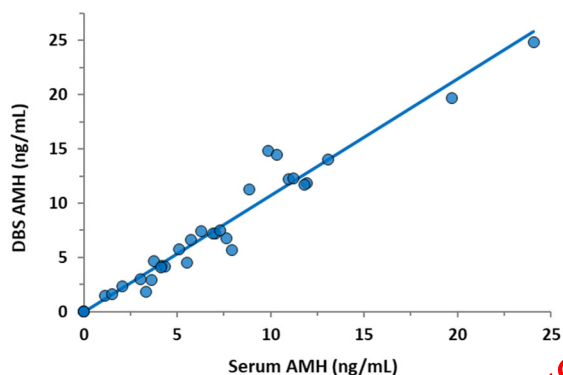

Proteolytic processing of anti-Mullerian hormone differs between human fetal testes and adult ovaries. *Molecular Human Reproduction* 2015; 21 (7): 571–582

2. di Clemente N, Jamin SP, Lugovskoy A, Carmillo P, Ehrenfels C, Picard JY, Whitty A, Josso N, Pepinsky RB, Cate RL. Processing of anti-mullerian hormone regulates receptor activation by a mechanism distinct from TGF-beta. *Mol Endocrinol.* 2010; 24 (11):2193-2206
3. Validation of Dried Blood Spot AMH ELISA: A Convenient Alternative to Venipuncture, Endocrine Society, 2016. Available: <https://endo.confex.com/endo/2016endo/webprogram/Paper26621.html>
4. HHS Publication, 5th ed., 2007. Biosafety in Microbiological and Biomedical Laboratories. Available <http://www.cdc.gov/biosafety/publications/bmbl5/BMBL5>
5. DHHS (NIOSH) Publication No. 78–127, August 1976. Current Intelligence Bulletin 13 - Explosive Azide Hazard. Available: <http://www.cdc.gov/niosh>.
6. Kricka L. Interferences in immunoassays – still a threat. *Clin Chem* 2000; 46: 1037–1038.

FOR RESEARCH USE ONLY

Not for use in diagnostic procedures.

The Ansh Labs logo is a trademark of Ansh Labs.

Manufactured by:
 Ansh Labs
 445 Medical Center Blvd.
 Webster, TX 77598-4217, U.S.A.

EXPECTED RESULTS

The expected results were determined by testing 20,738 specimens from a total of 13,555 healthy subjects whose blood was collected at multiple study sites in the US and Europe. Because AMH declines as ovarian reserve diminishes with age the values obtained using the AL-124-i ENSA are age stratified. AMH measurements are in pg/mL.

Age Range	Specimens Tested*	Mean	Mean SD	Lower Limit (2.5%)	Upper Limit (97.5%)
20-25	234	5018	331	1400	17356
26-30	542	4218	151	444	12774
31-35	985	3424	106	249	12213
36-40	1436	2375	67	36	8874
41-45	2723	881	21	<2	3866
46-50	5931	258	6	<2	1718
51-55	4546	46	2	<2	390
56-60	1972	6	3	<2	12
61-65	2369	2	<2	<2	3

* These 20,738 specimens were from a total of 13,555 individual women

Note: It is recommended that each laboratory should determine the reference range(s) for its own patient population. The results of this assay should be used in conjunction with other relevant and applicable clinical information.

REFERENCES

1. L.S. Mamsen, T.S. Petersen, J.V. Jeppesen, K. Mollgard, M.L. Grondahl, A. Larsen, E. Ernst, C. Oxvig6, A. Kumar, B. Kalra, and C.Y. Andersen.